TECHNICAL SERVICE BULLETIN 19-008



Date Issued: March 1, 2019

Subject: General MAPI Failure errors during MSDRx[®] Emailing Function

Explanation: This Technical Bulletin provides information concerning the emailing functionality of LGC Maine Standards' MSDRx[®] software and General MAPI Failure error messages when using Microsoft[®] Windows[®] Vista or newer operating systems.

MSDRX[®] emailing functionality:

The MSDRx[®] software includes functionality that allows the user to email data files and peer setup files generated by the software to LGC Maine Standards through Microsoft[®] Office Outlook[®] or other email programs.

General MAPI Failure error message:

When using the MSDRx[®] software, you may receive a "*General MAPI failure*" error message if Microsoft[®] Office Outlook[®] is open and the user clicks the Email button in the "E-mail data" section in MSDRx[®] after data files are selected. This error is caused by a Microsoft[®] security feature that was introduced with Windows[®] Vista operating system.

Resolution:

There are two options that will allow you to continue utilizing the MSDRx[®] software's emailing functionality if you receive the error mentioned above.

- 1. Close Outlook[®] before attempting to use the email functionality.
- 2. Open Outlook[®] with Administrator permissions before attempting to use the email functionality.

If these options fail, you can manually generate an email through Outlook[®] and attach the data files.

Please call LGC Maine Standards' Technical Support if you have additional questions regarding the emailing functionality of the MSDRx[®] software.

Contact Information:	Technical Support Department Maine Standards Company, LLC 221 US Route 1 Cumberland Foreside, ME 04110	
	Phone:	1-800-377-9684 ext. 1428
		+1-207-892-1300 ext. 1428
	Fax:	+1-207-892-2266
	E-mail:	msc.techsupport@LGCGroup.com